


Rtx[®]-CLPesticides Column

(30m, 0.25mm ID, 0.25µm-cat.# 11123)

Peak Identification

1. 2,4,5,6-tetrachloro-*m*-xylene
2. α -BHC
3. γ -BHC (lindane)
4. β -BHC
5. δ -BHC
6. heptachlor
7. aldrin
8. heptachlor epoxide
9. γ -chlordane
10. α -chlordane
11. 4,4'-DDE
12. endosulfan I
13. dieldrin
14. endrin
15. 4,4'-DDD
16. endosulfan II
17. 4,4'-DDT
18. endrin aldehyde
19. methoxychlor
20. endosulfan sulfate
21. endrin ketone
22. decachlorobiphenyl

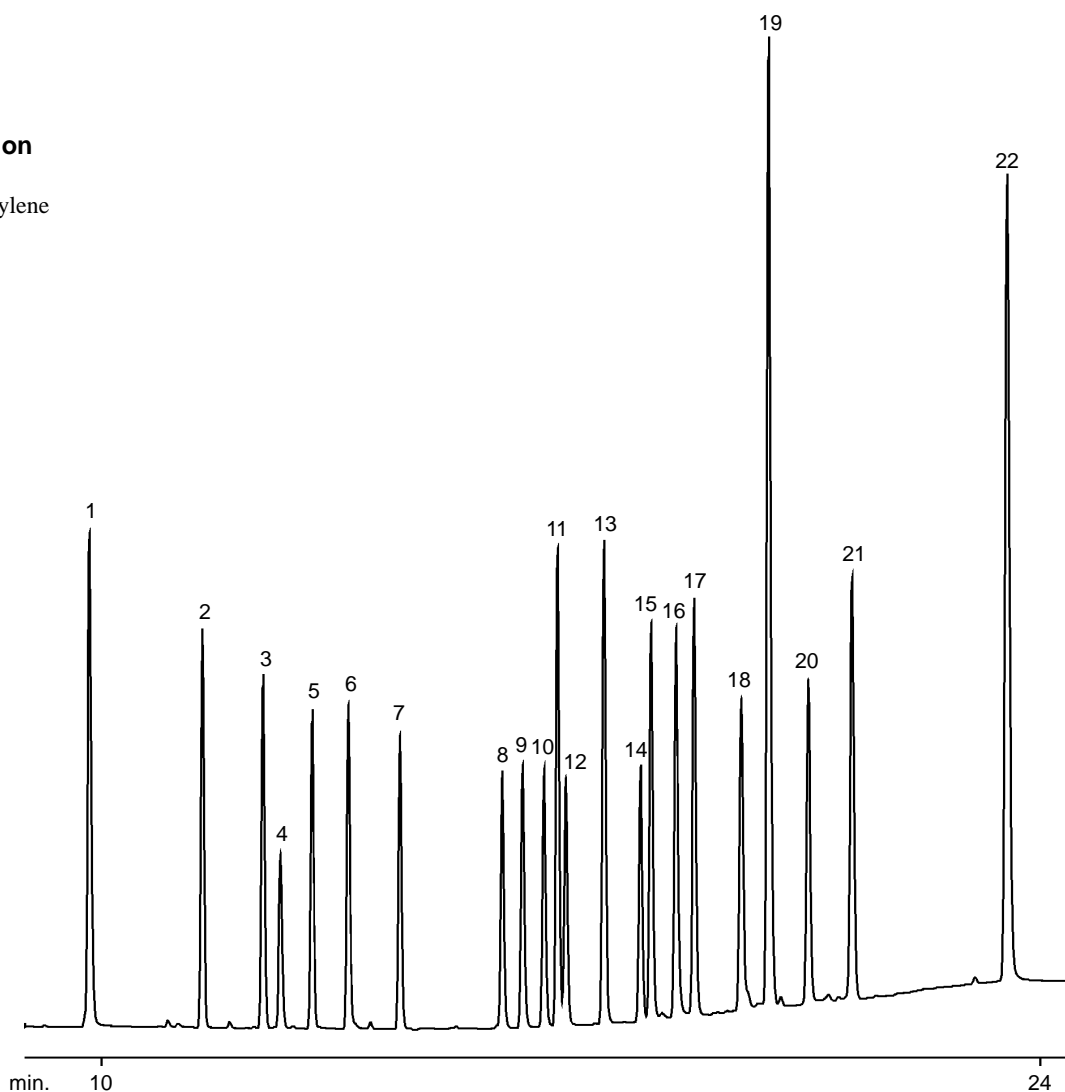

Run Conditions

On column concentration: 16-160 pg Organochlorine Pesticide Mix AB#2 (cat.# 32292); **Oven program:** 120°C (hold 1 min.) to 300°C @ 8.5°C/min. (hold 3 min.); **Injection port:** direct injection using a Uniliner[®] sleeve (cat.# 20335) at 200°C; **Detector:** ECD at 300°C with anode purge; **Dead time:** 2.4 minutes; **Head pressure:** 11.2 psi (constant); **Flow rate:** He @ 0.64mL/min. @ 120°C.

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



ISO 9001
#124-01 [002]
Rev. date: 5/01

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: dmannusrestek@compuserve.com