


# Rtx<sup>®</sup>-CLPesticides Column

(30m, 0.32mm ID, 0.50µm—cat.# 11139)

## Peak Identification

1. 2,4,5,6-tetrachloro-*m*-xylene
2. α-BHC
3. γ-BHC (lindane)
4. β-BHC
5. δ-BHC
6. heptachlor
7. aldrin
8. heptachlor epoxide
9. γ-chlordane
10. α-chlordane
11. 4,4'-DDE
12. endosulfan I
13. dieldrin
14. endrin
15. 4,4'-DDD
16. endosulfan II
17. 4,4'-DDT
18. endrin aldehyde
19. methoxychlor
20. endosulfan sulfate
21. endrin ketone
22. decachlorobiphenyl

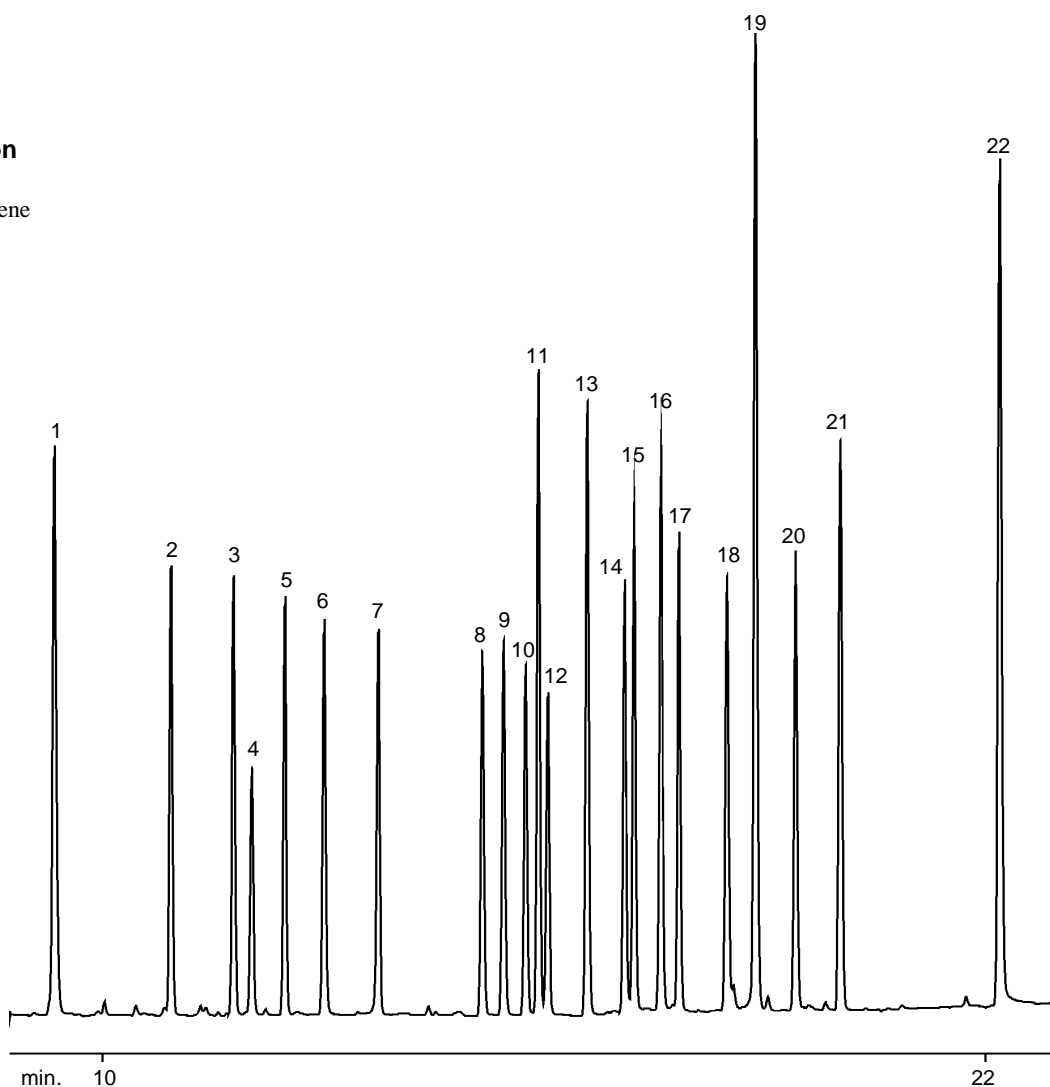

## Run Conditions

**On column concentration:** 16-160 pg Organochlorine Pesticide Mix AB#2 (cat.# 32292); **Oven program:** 120°C (hold 1 min.) to 300°C @ 9°C/min. (hold 10 min.); **Injection port:** direct injection using a Uniliner<sup>®</sup> sleeve (cat.# 20335) at 200°C; **Detector:** ECD at 300°C with anode purge; **Dead time:** 1.9 minutes; **Head pressure:** 8.7 psi (constant); **Flow rate:** He @ 1.3mL/min. @ 120°C.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.**



#124-01 [003]  
Rev. date: 5/01

**Restek USA:** 110 Benner Circle, Bellefonte, PA 16823  
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; [www.restekcorp.com](http://www.restekcorp.com)  
**Restek GmbH:** phone: (49) 06172 2797 0 • email: [RESTEK-GMBH@t-online.de](mailto:RESTEK-GMBH@t-online.de)  
**Restek France:** phone: (33) 01 60 78 32 10 • email: [restekfr@club-internet.fr](mailto:restekfr@club-internet.fr)  
**Thames Restek UK Ltd.:** phone: (44) 01753 624111 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)  
**Thames Restek Scotland Ltd.:** phone: 0870 241 1247 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)  
**Restek Ireland:** phone: (44)2 890 814 576 • email: [dmannusrestek@compuserve.com](mailto:dmannusrestek@compuserve.com)