

Inlet Liner Packing Tool

(cat.# 20339)

Overview

The inlet liner packing tool (Figure 1) is designed to pack the wool in the same position inside the liner to promote reproducibility from liner to liner. The tool also makes installation of the wool easy and hassle-free. The inlet liner packing tool is designed to be used with inlet liner IDs greater than 2mm.

Figure 1: The inlet liner packing tool.



Typical Packing Distances for Various Instruments*

Instrument	Packing Distance
Agilent 5890/6890/6850	27mm from open end
Varian 1177 1078/1079 1075/1077	37mm from open end 20mm from open end 34mm from open end
Carlo Erba 5000-6000 series 8000 series & Trace	35mm from open end 55mm from open end
Shimadzu 17A 14A	30mm from open end 35mm from open end
Perkin Elmer Autosys Sigma series Autosys XL	20mm from ground end 10mm from open end 50mm from open end
Finnigan 9001 and QCQ	27mm from open end

* Check the operating manual for packing distances for your particular instrument.

Instructions

1. Set the insertion distance on the inlet liner packing tool by loosening the knurled nut and sliding the scale/plunger until the proper distance is set for your application and instrument. After insertion distance is set, finger-tighten the knurled nut (do NOT overtighten).
2. Start the packing material into the liner (Figure 2).
3. Set the packing material to proper insertion distance by inserting the plunger into the liner until the base of the tool is flush with the liner end (Figure 3).

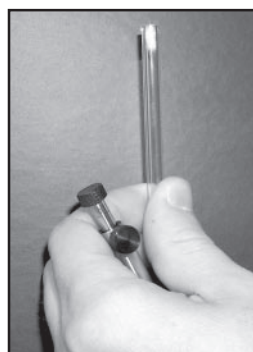


Figure 2: Start the packing material into the liner.



Figure 3: Set the packing material to proper distance using the liner packing tool.

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



ISO 9001
#200-00 [005]
Rev. date: 1/02/02

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: restekeurope@cs.com