


Instructions for Using Alumaseal™ Ferrules

Cat.# 21472, 21473, 21474

1. Place column nut onto capillary tubing.
2. Place back of Alumaseal™ ferrule onto capillary tubing as shown in figure.
3. Place front of Alumaseal™ ferrule onto capillary tubing.
4. Tighten nut until capillary tubing cannot be moved.

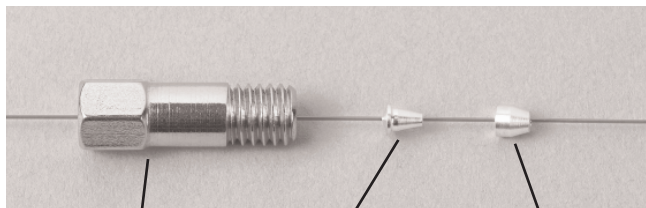

Column Nut

Back of ferrule

Front of ferrule



110 Benner Circle
Bellefonte, PA 16823-8812
814-353-1300 H 800-356-1688
fax: 814-353-1309
www.restekcorp.com

#201-05 [001]
©2004 Restek Corporation