


Moisture Control System (MCS) By-Pass Line

For use with Tekmar 3000 Purge & Trap

(Includes 1/16" OD x 0.020" ID Silcosteel® tubing; two 1/16" stainless steel Valco® nuts; and two stainless steel Valco® ferrules—cat.# 21035)

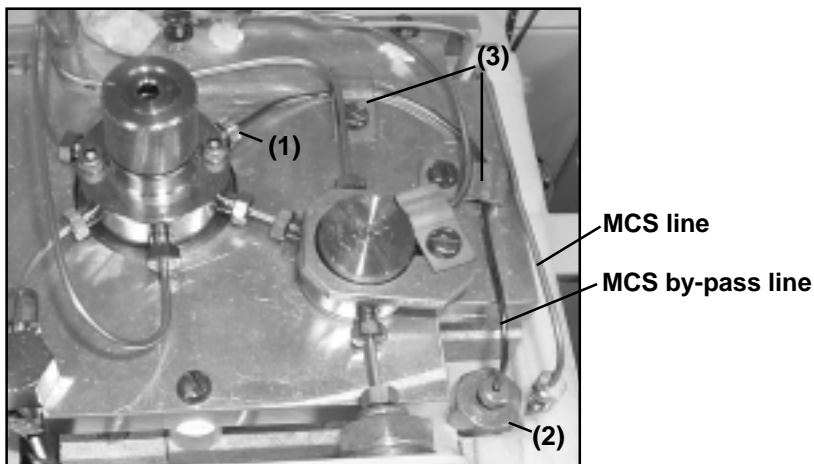

1. Remove existing MCS by-pass line from the instrument by removing the two 1/16-inch nuts from the instrument (1 and 2).
2. Place one nut and one ferrule onto one end of the MSC by-pass line and install onto (1).
3. Run the MCS by-pass line through the existing clamps (3) and tighten.
4. Place one nut and one ferrule onto the other end of the MCS by-pass line and install onto (2).

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



#203-02 [022]
Rev. date: 6/16

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: dmannusrestek@compuserve.com