

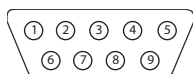


Remote Start Cable

Agilent 5890 GC to Agilent 35900C Interface
 Agilent 5890 GC to Agilent 7673A Autosampler
 (cat.#s 20660, 20661—replaces Agilent #35900-60700)

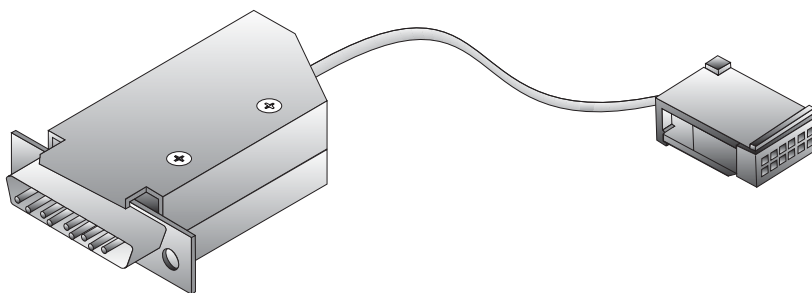

Connector 1
front view

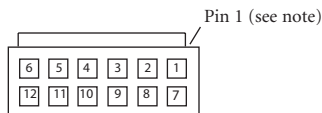

Connector 2
front view

Conn. 1 Pin	Conn. 2 Pin	Color	Function
1	11	Black	Ground
3	9	Red	Start sampler or interface
7	4	White	Ready
	1, 3		Jumper
	6, 8		Jumper
	10, 12		Jumper

Note: Connector 2 pin numbers correspond to the pin numbers on an Agilent 5890 GC. Pin 1 is stamped Pin 1 on connector.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4
 (or your Restek representative) if you have any questions about this
 product or any other Restek product.**



#203-03 [006]
 Rev. date: 01/07

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823
 phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • www.restek.com
Restek France • phone: 33 (0)1 60 78 32 10 • e-mail: restek@restekfrance.fr
Restek Ireland • phone: 44 2890 814576 • e-mail: resteurope@aol.com
Thames Restek U.K. LTD • phone: 44 1494 563377 • e-mail: sales@thamesrestek.co.uk
Restek GmbH • phone: +49 (0) 6172 2797 0 • e-mail: info@restekgmbh.de