


# Remote Start Cable

Agilent 3396 Integrator to Remote Start  
(cat.#s 20655, 20656—replaces Agilent #03394-60540)

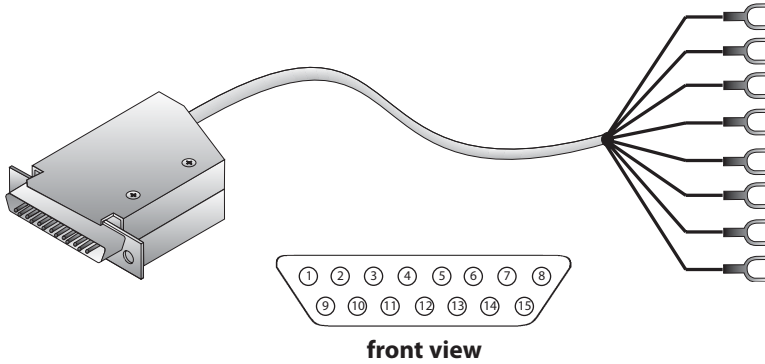

Name	Pin	Color	Function
Start in	3	Green	Start integrator
Stop in	6	White	Stop integrator
Start out 1	1	Black	Start external device
Start out 2	13	Orange	Start external device
Ready in	5	Red	Indicate readiness to integrator
Ready out	14	Blue	Indicate integrator readiness
Ground	15	Brown	Ground

## Description

- Start in:** Starts integrator when momentarily grounded.  
**Stop in:** Stops integrator when momentarily grounded.  
**Start out:** One-second relay contact closure between Pins 1 and 13.  
**Ready out:** TTL high (5 volt) when integrator ready light is lit.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4  
(or your Restek representative) if you have any questions about this  
product or any other Restek product.**



#203-03 [010]  
Rev. date: 01/07

**Restek U.S.** • 110 Benner Circle • Bellefonte, PA 16823

phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • [www.restek.com](http://www.restek.com)

**Restek France** • phone: 33 (0)1 60 78 32 10 • e-mail: [restek@restekfrance.fr](mailto:restek@restekfrance.fr)

**Restek Ireland** • phone: 44 2890 814576 • e-mail: [resteurope@aol.com](mailto:resteurope@aol.com)

**Thames Restek U.K. LTD** • phone: 44 1494 563377 • e-mail: [sales@thamesrestek.co.uk](mailto:sales@thamesrestek.co.uk)

**Restek GmbH** • phone: +49 (0) 6172 2797 0 • e-mail: [info@restekgmbh.de](mailto:info@restekgmbh.de)