



Optional Split Vent Trap Assembly

(Kit-cat.# 23031; Replacement Trap-cat.# 23032)


Installation Instructions: For use with Agilent 6890/6850 GC's.

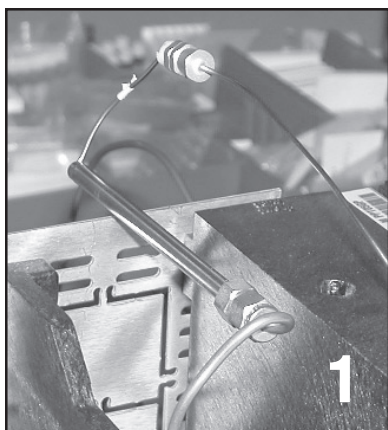


Remove top cover from the rear of the GC. Locate the pencil trap inside the GC (the trap is located at the rear of the GC on the injector side)

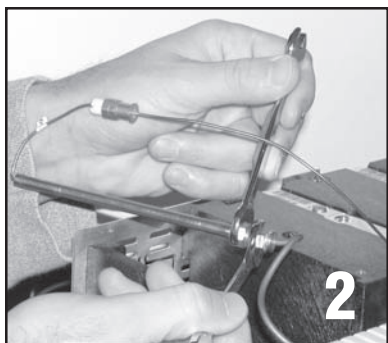


Remove the trap by loosening and removing the 1/8" nut.

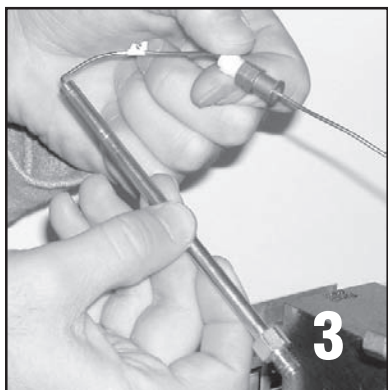


Loosen and remove the finger tight nut from the trap.

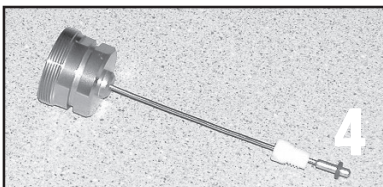


Place the new plastic finger tight nut and o-ring onto the new optional split vent trap.

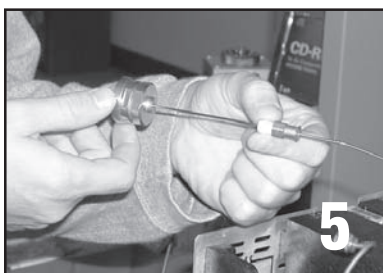


Tighten the new finger tight nut and o-ring assembly into the existing GC connection.

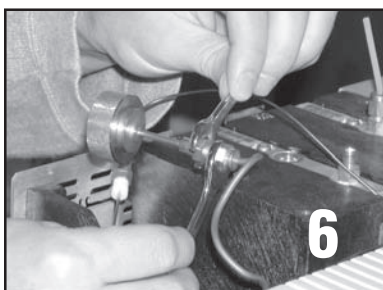


Tighten the other half of the new optional split trap assembly onto the existing 1/8" fitting. (Restek recommends that you replace the 1/8" nut and ferrule in order to ensure a leak free connection).

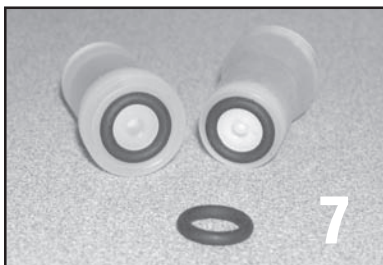

Place the o-rings onto the ends of the trap.

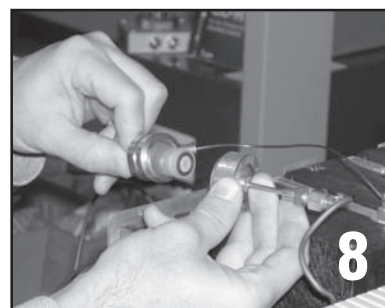

Place the trap into the new optional split vent trap housing.

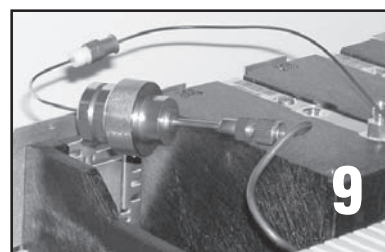

Tighten the two fittings together by tightening the knurled nut. (Do not use wrenches or other instruments to tighten the fitting)

**Call Technical Service at
800-356-1688 or
814-353-1300, ext. 4
(or your local Restek
representative)
if you have any questions
about this product or any
other Restek product.**

RESTEK
Turning Visions Into Reality

ISO 9001
#20303 [022]
Rev. date: 1/03

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01494 563377 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: dmannusrestek@compuserve.com