



# Capillary Inlet Adaptor Fitting for Agilent GCs

cat.# 27184

## Installation Instructions

(To be used with single hole ferrules only)

1. Cool GC and remove the old inlet fitting.
2. Assemble as shown below and screw tightly into the GC.
3. Insert the column 30-32mm from the bottom of the  $\frac{1}{16}$ -inch nut.
4. Leak check with a Restek Leak Detector (cat.# 22839).

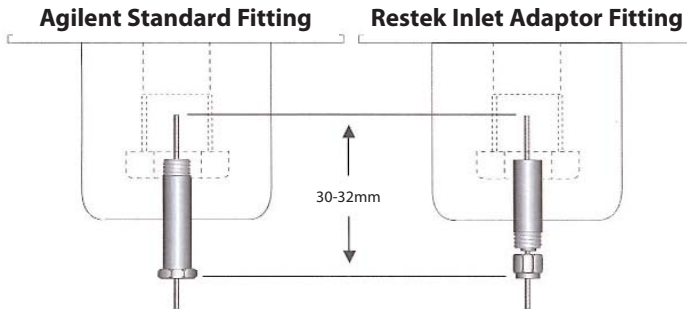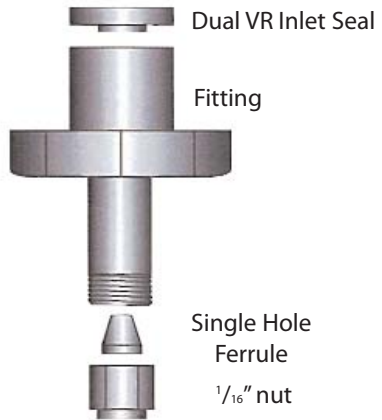

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.



Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823

phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • [www.restek.com](http://www.restek.com)

Turning Visions Into Reality

Restek France • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: [restek@restekfrance.fr](mailto:restek@restekfrance.fr)

Restek Ireland • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: [restekeurope@aol.com](mailto:restekeurope@aol.com)



Thames Restek U.K. LTD • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: [sales@thamesrestek.co.uk](mailto:sales@thamesrestek.co.uk)

#203-03 [037]

Restek GmbH • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: [info@restekgmbh.de](mailto:info@restekgmbh.de)

Rev. date: 1/09