


ECD Vent Trap Kit

(Includes ECD Vent Trap and 9-inch piece of copper tubing—cat.# 22017)

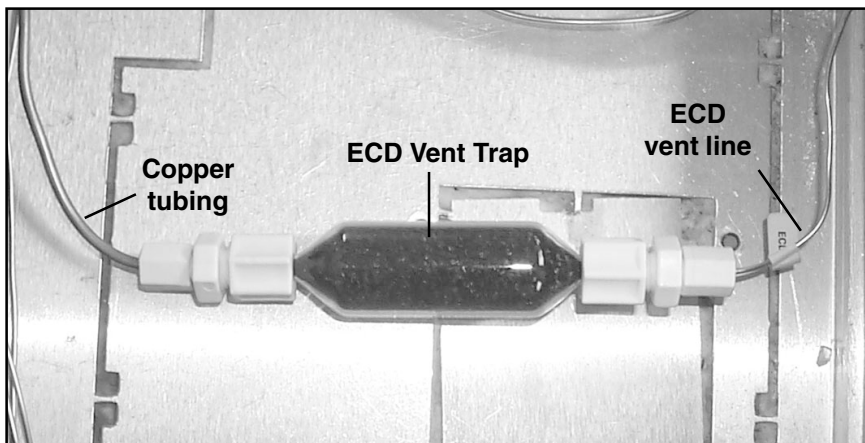

1. Connect the 9-inch piece of copper tubing to one end of the ECD vent trap.
2. Connect the other end of the ECD vent trap to the existing ECD vent line.
3. Position the ECD vent trap in the GC as desired. Avoid placing the trap in locations where it could be broken.

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



#204-08 [009]
Rev. date: 1/01

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: dmannusrestek@compuserve.com