

Outlet Cap & Gold Seal Assembly Tool

for Agilent 1100 HPLC Systems



**EASILY INSTALL
GOLD SEAL INTO
OUTLET CAP**



Put Outlet Cap on male part of Assembly Tool.



Push the Assembly Tool together.



Hold onto Outlet Cap and pull Assembly Tool apart.



Put Gold Seal over pin on male part of Assembly Tool.



Push the Assembly Tool together and press the Gold Seal in the Outlet Cap.



Pull the Assembly Tool apart and remove assembled Outlet Cap and Gold Seal.

Description

Outlet Cap and Gold Seal Assembly Tool for Agilent 1100 HPLC Systems

qty.

ea.

cat.#

24989

Restek Replacement Parts for Agilent HPLC Systems

• Meet or exceed OEM performance.

Description

Description	Model #	Similar to Agilent part #	qty.	cat.#
Seal, Gold Disk (outlet)	1050 & 1100	5001-3707	ea.	25467
Outlet Cap	1050 & 1100	5062-2485	4-pk.	25139
Outlet Cap & Gold Seal Assembly	1050 & 1100	—	2-pk.	25140



Chromatography Products

Lit. Cat.# 580159

© 2007 Restek Corporation.

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823 • 814-353-1300 • 800-356-1688 • fax: 814-353-1309 • www.restek.com

Restek France • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: restek@restekfrance.fr

Restek Ireland • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: restekeurope@aol.com

Thames Restek U.K. LTD • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: sales@thamesrestek.co.uk

Restek GmbH • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: info@restekgmbh.de

