



# Trident™ In-Line 2cm Guard Cartridge Holder

(cat.#s 25060 and 25061)

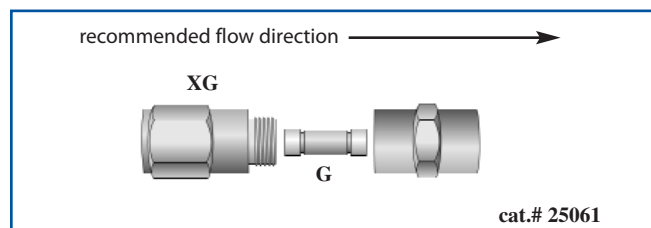

The **Trident™ in-line guard cartridge holder** can be used with almost all HPLC columns by connecting it with a short piece of 1/16" tubing, appropriate nuts and ferrules, or finger-tight fittings. To use, remove the XG fitting, insert the appropriate guard cartridge (G), and then reinstall the XG fitting. When tightening the fitting, turn it no more than 1/8-turn past finger-tight, as there is no need to overtighten this connection.

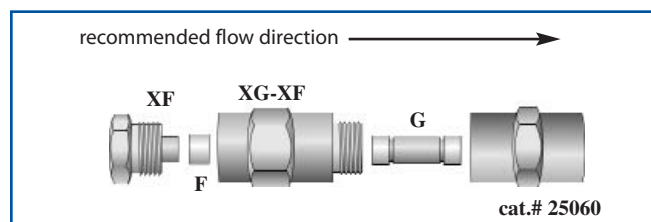

The **Trident™ in-line guard cartridge holder with filter** can be used with almost all HPLC columns by connecting it with a short piece of 1/16" tubing, appropriate nuts and ferrules, or fingertight fittings. To use, remove the XG-XF fitting, insert the appropriate guard cartridge (G), and then reinstall the XG-XF fitting. When tightening the fitting, turn it no more than 1/8-turn past finger-tight, as there is no need to overtighten this connection. To replace the cap frit (F), remove the XF fitting, pull off the old frit, push a new frit into place, and then reinstall the XF fitting. When tightening the XF fitting, turn no more than 1/8-turn past finger-tight.

Restek Trademarks: Trident, Restek logo, Turning Visions Into Reality  
Other Trademarks: PEEK (Victrex plc).

**For a complete listing of guard cartridges, please visit [www.restek.com](http://www.restek.com).**

## Trident™ In-line 2cm Guard Cartridge Holders

- 2cm Guard Cartridge Holder: cat.# 25061
- 2cm Guard Cartridge Holder with Filter: cat.# 25060

## Replacement parts for the Trident™ In-line Column Protection System

- 1/16" OD PEEK™ Tubing:
  - 0.0025" ID Natural cat.# 25320
  - 0.005" ID Red Stripe cat.# 25065
  - 0.007" ID Yellow Stripe cat.# 25066
  - 0.010" ID Blue Stripe cat.# 25067
  - 0.020" ID Orange Stripe cat.# 25068
- PEEK™ Column Connector: cat.# 25015
- PEEK™ Column Plug: cat.# 25016
- XF Fitting: cat.# 25024
- XG Fitting 2cm: cat.# 25063
- XG-XF Fitting 2cm: cat.# 25062
- Cap Frits 4mm, 2.0µm: cat.# 25022, 5-pk.
- Cap Frits 4mm, 0.5µm: cat.# 25023, 5-pk.
- Cap Frits 2mm, 2.0µm: cat.# 25057, 5-pk.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.**



ISO 9001:2000  
cert. # FM80397 #730-10-002  
Rev. date: 05/07

**Restek U.S.** • 110 Benner Circle • Bellefonte, PA 16823  
phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • [www.restek.com](http://www.restek.com)  
**Restek France** • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: [restek@restekfrance.fr](mailto:restek@restekfrance.fr)  
**Restek Ireland** • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: [restekeurope@aol.com](mailto:restekeurope@aol.com)  
**Thames Restek U.K. LTD** • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: [sales@thamesrestek.co.uk](mailto:sales@thamesrestek.co.uk)  
**Restek GmbH** • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: [info@restekgmbh.de](mailto:info@restekgmbh.de)