



# Trident™ Direct In-Line Filter and 2cm Guard Cartridge Holders

(cat.#'s 25085 and 25086)

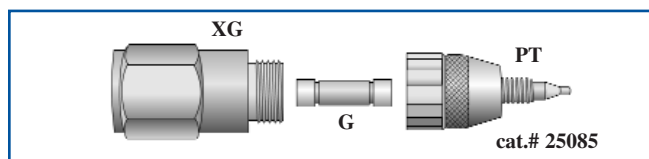

The **Trident™ Direct in-line guard cartridge holder** can be used with all HPLC columns that have CPI-style\* end fittings. To use, remove the XG fitting, insert the appropriate guard cartridge (G), and then reinstall the XG fitting. When tightening the fitting, turn it no more than 1/8-turn past finger-tight, as there is no need to overtighten this connection. Then simply insert the PEEK™ tip (PT) into the column inlet and finger tighten.

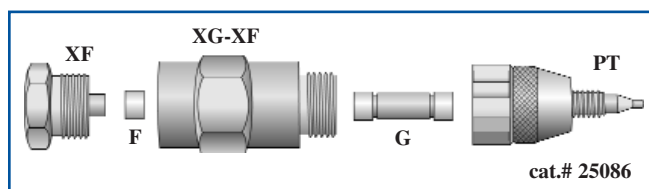

The **Trident™ Direct in-line guard cartridge holder with filter** can be used with all HPLC columns that have CPI-style\* end fittings. To use, remove the XG-XF fitting, insert the appropriate guard cartridge (G), and reinstall the XG-XF fitting. When tightening the fitting, turn it no more than 1/8-turn past finger-tight, as there is no need to overtighten this connection. Then simply insert the PEEK™ tip (PT) into the column inlet and finger tighten. To replace the cap frit (F), remove the XF fitting, pull off the old frit, push a new frit into place, and then reinstall the XF fitting. When tightening the XF fitting, turn no more than 1/8-turn past finger-tight.

**For a complete listing of guard cartridges, please visit [www.restek.com](http://www.restek.com).**

## Trident™ Direct High-Pressure Filter and Guard Cartridge Holders

- 2cm Guard Cartridge Holder: cat.# 25085
- 2cm Guard Cartridge Holder with Filter: cat.# 25086

## Replacement parts for the Trident™ Direct Column Protection System

- PEEK™ Tip for CPI Fittings: cat.# 25087
- PEEK™ Tip for Waters Fittings: cat.# 25088
- XF Fitting: cat.# 25024
- XG Fitting 2cm: cat.# 25063
- XG-XF Fitting 2cm: cat.# 25062
- Cap Frits 4mm, 2.0µm: cat.# 25022, 5-pk.
- Cap Frits 4mm, 0.5µm: cat.# 25023, 5-pk.
- Cap Frits 2mm, 2.0µm: cat.# 25057, 5-pk.

\* The standard PEEK™ tip that comes with the Trident™ Direct system is compatible with Parker®, Swagelok®, Upchurch®, Valco®, and other CPI-style fittings. To use the Trident™ Direct system with a Waters® style fitting, the tip must be replaced with cat.# 25088.

Restek Trademarks: Trident, Restek logo, Turning Visions Into Reality

Other Trademarks: PEEK (Victrex plc), Parker (Parker Instrumentation Division), Swagelok (Swagelok Company), Upchurch (Upchurch Scientific), Valco (Valco Instruments Co, Inc.), Waters (Waters Associates, Inc.).

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.**



#730-10-004  
Rev. date: 05/07

**Restek U.S.** • 110 Benner Circle • Bellefonte, PA 16823  
phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • [www.restek.com](http://www.restek.com)

**Restek France** • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: [restek@restekfrance.fr](mailto:restek@restekfrance.fr)

**Restek Ireland** • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: [restekeurope@aol.com](mailto:restekeurope@aol.com)

**Thames Restek U.K. LTD** • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: [sales@thamesrestek.co.uk](mailto:sales@thamesrestek.co.uk)

**Restek GmBH** • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: [info@restekgmbh.de](mailto:info@restekgmbh.de)