


LC Cartridge Holders


(cat.# 25298, 25299)

For Direct Connections

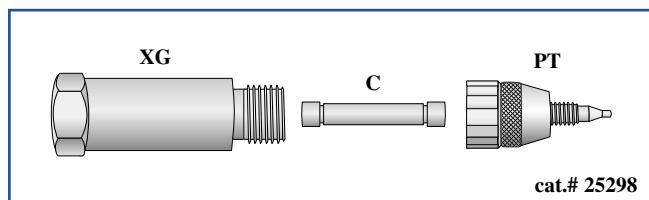

The Restek Direct Connect Fast LC Cartridge Holder with filter can be used with any instrument that has CPI-style* fittings. To use, remove the XG fitting, insert the fast LC cartridge (C), and reinstall the XG fitting. Tighten the fitting no more than $\frac{1}{8}$ -turn past finger-tight, as there is no need to over-tighten this connection. Then simply insert the PEEK® tip (PT) into the detector, or connect to capillary tubing, using a union, and finger tighten.

For In-Line Connections

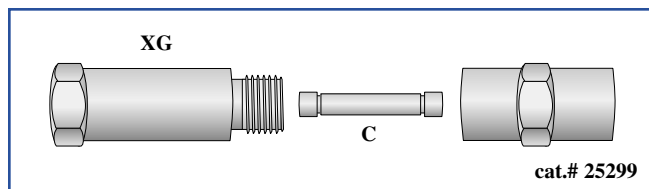

To use, remove the XG fitting, insert the fast LC cartridge (C), and reinstall the XG fitting. Tighten the fitting no more than $\frac{1}{8}$ -turn past finger-tight, as there is no need to over-tighten this connection. Then simply connect the holder to the instrument like a conventional column.

Fast LC Cartridges

Phase	Particle Size	Dimensions	cat.#
Ultra Cyano	3µm	30 x 2.1mm	91063320
Ultra Cyano	3µm	30 x 4.0mm	91063340
Ultra C18	3µm	30 x 2.1mm	91743320
Ultra C18	3µm	30 x 4.0mm	91743340
Ultra PFP	3µm	30 x 2.1mm	91763320
Ultra PFP	3µm	30 x 4.0mm	91763340
Ultra Aqueous C18	3µm	30 x 2.1mm	91783320
Ultra Aqueous C18	3µm	30 x 4.0mm	91783340

Fast LC Cartridge Holders

- for Direct Connections: cat.# 25298
- for In-Line Connections: cat.# 25299

Replacement Parts for the Direct Connect Fast LC System

- PEEK® Tip for CPI Fittings: cat.# 25087
- PEEK® Tip for Waters Fittings: cat.# 25088

*The standard PEEK® tip included with the Restek Fast LC System is compatible with Parker®, Swagelok®, Upchurch®, Valco®, and other CPI-style fittings. To use the Restek Fast LC System with a Waters® style fitting, the tip must be replaced with cat.# 25088.

For trademark attributions please refer to our current catalog.

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



ISO 9001
#730-10 [005]
Rev. date: 1/31/03

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • fax: (49) 06172 2797 77
Restek France: phone: 01 60 78 32 10 • fax: 01 60 78 70 90
Thames Restek UK Ltd.: phone: 01494 563377 • fax: 01494 564990
Restek Ireland: phone: 44 28 9081 4576 • fax: 44 28 9081 4576