

2-in-1 Filter/O-Ring Insertion Tool Kit

for ASE® 100/200/300 Extraction Cells

Inserting a filter (ASE® 100/200/300 systems), using Filter Insertion Attachments on Resprep™ Tool Handle.



Screw the appropriate size attachment onto the end of the Resprep™ Tool Handle.



Place a filter at the top of the extraction cell.



Push the filter to the bottom of the extraction cell.



2-in-1 Filter/O-Ring Insertion Tool Kit

Inserting an O-Ring (ASE® 100/200/300 systems), using Resprep™ Tool Handle.



Place the O-ring over the tip of the tool.



Insert the tool into the center hole of the extraction cell cap.



Press the tool firmly inside the cap until the O-ring snaps into place.



Resprep™ Tool Handle



Filter Insertion Attachments

Description

Description	qty.	cat.#
2-in-1 Filter/O-Ring Insertion Tool Kit for ASE® 100/200/300 (includes Resprep™ Tool Handle and Filter Insertion Attachments)	kit	26181
Resprep™ Tool Handle for ASE® 100/200/300	ea.	26182
Filter Insertion Attachments for ASE® 100/200/300 (1mL, 5mL, 11mL, 33mL)	4-piece set	26183

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.



#803-04-001
Rev. date: 10/07

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restek.com
Restek GmbH: phone: (49) 06172 2797 0 • fax: (49) 06172 2797 77
Restek France: phone: 01 60 78 32 10 • fax: 01 60 78 70 90
Thames Restek UK Ltd.: phone: 01494 563377 • fax: 01494 564990
Restek Ireland: phone: 44 2890 814576 • fax: 44 2890 814576