

Tubing Assembly for ASE® 200 Systems

cat.# 26251

Restek's Tubing Assembly for ASE® 200 Systems eliminates the need for an adaptor fitting and allows 1/32" tubing to be inserted directly into either static or pressure relief valves. This simplifies installation and removes a potential source of leaks.

Installation Instructions

1. Remove the old tubing assembly and adaptor from the static or pressure relief valve. The adaptor is not needed when using Restek's tubing assembly for ASE® 200 Systems (Figure 1).
2. Install the 1/16" tubing side of Restek's Tubing Assembly for ASE® 200 Systems directly into the static or pressure relief valve, whichever applies (Figure 2).
3. Install the 1/32" tubing side of Restek's Tubing Assembly for ASE® 200 Systems into the upper or lower AutoSeal tip, whichever applies.

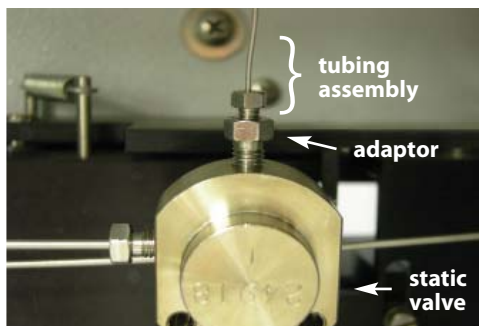


Figure 1

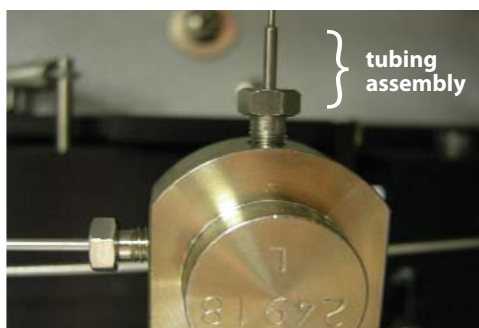


Figure 2

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.



Chromatography Products
© 2009 Restek Corporation



#803-04 [002]
Rev. date: 8/09

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823
phone: 814-353-1300 • 800-356-1688 • fax: 814-353-1309 • www.restek.com
Restek France • phone: +33 (0)1 60 78 32 10 • e-mail: restek@restekfrance.fr
Restek GmbH • phone: +49 (0)6172 2797 0 • e-mail: info@restekgmbh.de
Restek Ireland • phone: +44 (0)2890 814576 • e-mail: restekeuropa@aol.com
Restek Japan • phone: +81 (3)6459 0025 • e-mail: ryosei.kanaguchi@restek.com
Thames Restek U.K. LTD • phone: +44 (0)1494 563377 • e-mail: sales@thamesrestek.co.uk