

# Capillary Column Caps

(cat.# 21044)

---

## When to use Capillary Column Caps

- Anytime you want to store your capillary column.

## Why use Capillary Column Caps

- To replace using septa and avoiding septa particles from getting into your column.
- To create an airtight seal, which eliminates air and moisture from entering the column.

## How to use Capillary Column Caps

- Remove the capillary column from the GC.
- Make sure the column has square cuts on both ends.
- Insert the injection port end into the blue cap.
- Insert the detector port end into the clear cap.

The column is now ready for storage.

***Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.***

---



Turning Visions Into Reality



#202-04 [002]  
Rev. date: 5/01

**Restek USA:** 110 Benner Circle, Bellefonte, PA 16823

phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; [www.restekcorp.com](http://www.restekcorp.com)

**Restek GmbH:** phone: (49) 06172 2797 0 • email: [RESTEK-GMBH@t-online.de](mailto:RESTEK-GMBH@t-online.de)

**Restek France:** phone: (33) 01 60 78 32 10 • email: [restekfr@club-internet.fr](mailto:restekfr@club-internet.fr)

**Thames Restek UK Ltd.:** phone: (44) 01753 624111 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)

**Thames Restek Scotland Ltd.:** phone: 0870 241 1247 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)

**Restek Ireland:** phone: (44)2 890 814 576 • email: [dmannusrestek@compuserve.com](mailto:dmannusrestek@compuserve.com)