


FID Capillary Column Adapter for PerkinElmer Autosys™ XL

For use with 1/16" compression-style nuts, cat.# 22609

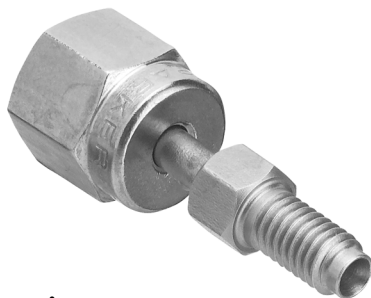

Installation Instructions

1. Cool the GC FID and remove the FID fitting.
2. Install the FID Capillary Column Adapter.
3. Install the 1/16" nut onto column, followed by a 1/16" capillary column ferrule.

Note: Always point the end of the column downward when placing a ferrule onto a column, to prevent any ferrule shards from entering the column.

4. Cut the column, using a scoring wafer. Verify that you have a square cut on the end of the column.
5. Insert the column 67-69 mm from the bottom of the 1/16" nut and tighten the nut ¼ turn past finger tight.
6. Check the fittings for leaks, using an electronic leak detector.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4
(or your Restek representative) if you have any questions about this
product or any other Restek product.**



#203-02 [027]
Rev. date: 5/05

Restek USA: 110 Benner Circle, Bellefonte, PA 16823

phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restek.com

Restek GmbH: phone: +49 (0) 6172 2797 0 • email: info@restekgmbh.de

Restek France: phone: 33 (0)1 60 78 32 10 • email: restekfr@club-internet.fr

Thames Restek UK Ltd.: phone: 44 1494 563377 • email: Sales@Thamesrestek.co.uk

Restek Ireland: phone: 44 2890 814576 • email: restekeuropa@aol.com