

Hot Swap Column Nut for Agilent GCs

For use with standard 1/16-inch ferrules

cat.# 22347



Recommended for use with graphite ferrules.

Installation instructions:

1. Verify the GC oven temperature is 50°C or less.
(Note, the Hot Swap Capillary Column Nut cools with the GC oven temperature. If the oven door is left open with the fan off for longer than 1 minute, the Hot Swap Capillary Column Nut will begin to heat to the inlet set temperature.)
2. Insert the capillary column through the Hot Swap Capillary Column Nut and ferrule.
3. Adjust the installation distance according to the manufacturer's recommendation.
4. Thread the Hot Swap Capillary Column Nut onto the injection port and turn until finger tight.
5. Leak check with a Restek Leak Detector (cat.# 22839).

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.



Turning Visions Into Reality



Rev. date: 12/08

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823

phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • www.restek.com

Restek France • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: restek@restekfrance.fr

Restek Ireland • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: restekeurope@aol.com

Thames Restek U.K. LTD • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: sales@thamesrestek.co.uk

Restek GmbH • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: info@restekgmbh.de